


## VenturiOne™ Trial Activation Process

There is a simple two step process to follow in order to activate the VenturiOne™ trial software;

1. We need to obtain information from the PC that you intend to run the software on, this is in the form of a license request file
2. Upon receipt of this information we will send by return an activation key code that will permit the software to run

Please see instructions below:

Upon installation of the software a licensing dialog box will appear as shown below.

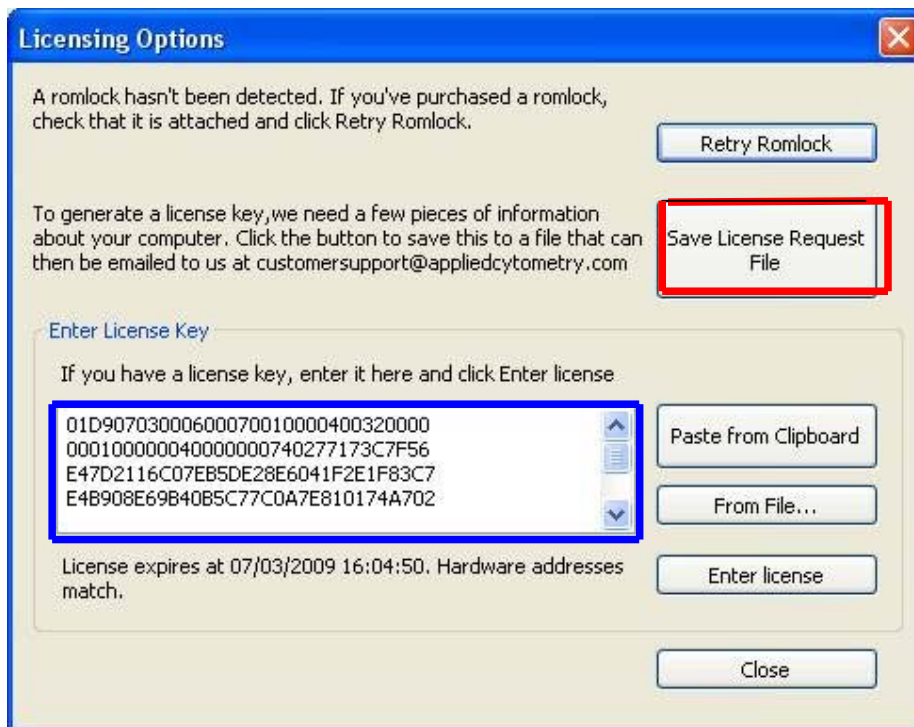

- As prompted, save the license request file, indicated by the red box (We recommend that you save this to your desk top for easy access)
- Email this file as an attachment to [customersupport@appliedcytometry.com](mailto:customersupport@appliedcytometry.com)
- By return a licence key will be emailed back to you
- When you receive the license key paste the code into the appropriate location in the licensing option dialog box, indicated by the blue box
- Click on 'Enter license' and the VenturiOne™ software should now be activated

If you need any further assistance, please do not hesitate to contact us at [customersupport@appliedcytometry.com](mailto:customersupport@appliedcytometry.com)